


## No. 4

### Schlieren Imaging of Shock Wave from Trigger Air Spark Gap

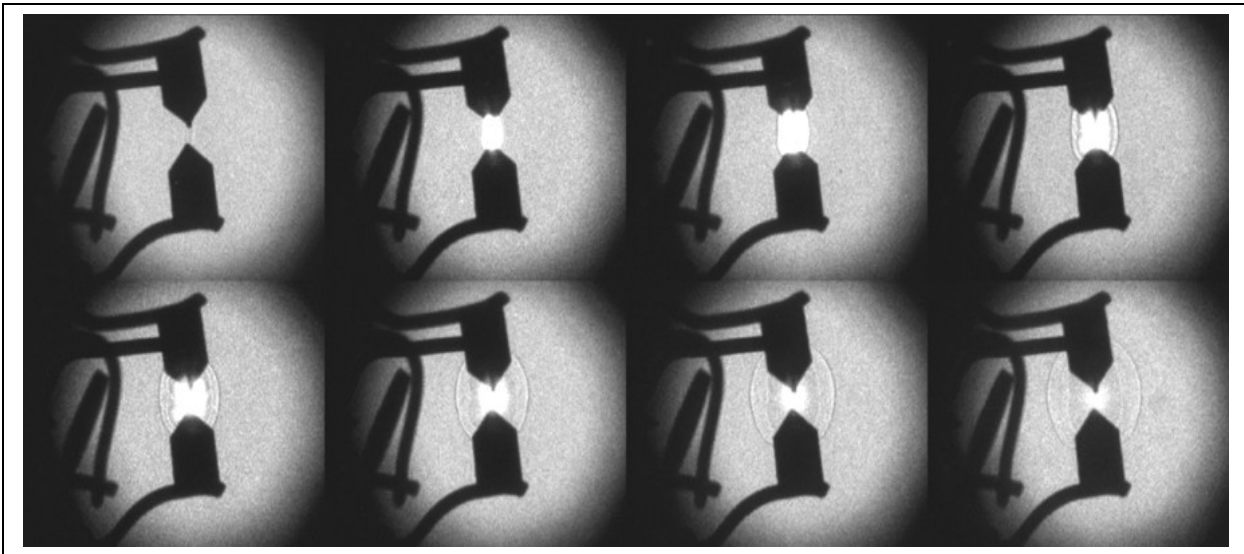

#### IMAGING PARAMETERS

SIM8 Camera running at 5 Million fps, 35 ns  
Exposures with Pseudo Schlieren Optical system.  
The field of view is about 10cm x 8cm.

#### OVERVIEW OF EXPERIMENT

The Air Spark Gap with 0.75 uF Capacitor  
and 1.5KV Charge Voltage across the 2 mm air gap.  
Two of the SIM's output triggers were used to  
synchronize the triggering of the spark gap and to  
pre-trigger the flash lamp used for the Schlieren  
light source.

#### EQUIPMENT PARAMETERS

A 1.5 msec duration xenon flash lamp directly coupled to a 0.125 inch diameter fibre light guide which was the point source. A 200 mm f2.5 lens was used to collimate the point source into a parallel beam, The Objective lens, a Nikon 85 mm f1.8 with extension tubes, was used to both focus the event into the camera at about 1:1 magnification. With the lens position away from the camera using the extension tube, the parallel light focuses at the infinity focus point of the objective lens, which then re-diverges to produce the background light on the image plane to produce the Schlieren imaging effect.

#### UK (Head Office / Factory)

6 Harvington Park, Pitstone Green  
Business Park, Pitstone.  
LU7 9GX England

**+44 (0) 1442 827728**

#### USA

Specialised Imaging Inc.  
40935 County Center Dr. Suite D  
Temecula, CA 92591, USA

**+1 951-296-6406**

#### GERMANY

Hauptstr. 10,  
82275 Emmering  
Germany

**+49 8141 666 89 50**



FM 87429 ISO9001:2015

